

## Lipofectamine® 2000 Reagent

Cat. nos. 11668-030	Size: 0.3 mL	Store at 4°C (do not freeze)
11668-027	0.75 mL	
11668-019	1.5 mL	
11668-500	15 mL	
Pub. Part No. 11668.2k.pps	Pub. No. MAN0000995	

### Description

- Lipofectamine® 2000 Transfection Reagent is a proprietary formulation for transfecting nucleic acids (DNA, RNA, and mRNA) into a wide range of eukaryotic cells.
- Nucleic acid-Lipofectamine® 2000 complexes can be added directly to cells in culture medium, in the presence or absence of serum/antibiotic.
- It is not necessary to remove complexes or change/add medium after transfection.

### Important Guidelines for Transfection

- Use Opti-MEM® I Reduced Serum Medium (Cat. no. 31985-062) to dilute Lipofectamine® 2000 Transfection Reagent and nucleic acids.
- The amount of Lipofectamine® 2000 Reagent required for successful transfection varies depending on the cell type and passage number. Start any new transfection by testing the recommended four concentrations of Lipofectamine® 2000 Reagent to determine an optimum amount.
- Prepare high-quality plasmid DNA at 0.5–5 µg/µL in deionized water or TE buffer. A GFP (green fluorescent protein) plasmid can be used to determine transfection efficiency.
- For additional information, and protocols on transfecting suspension cells refer to the manual at [www.lifetechnologies.com/transfection](http://www.lifetechnologies.com/transfection).

Day 0


Cells



Lipofectamine® 2000 Reagent



DNA

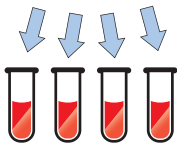

Day 1

Day 2-4

Seed cells so they are 70-90% confluent at the time of transfection.

Dilute 4 amounts of Lipofectamine® 2000 Reagent in Opti-MEM® Medium.

Dilute DNA in Opti-MEM® Medium.

Add diluted DNA to each tube of diluted Lipofectamine® 2000 Reagent (1:1 ratio).

Incubate.

Add DNA-reagent complex to cells.

Visualize/analyze transfected cells.

## Plasmid Transfection

Transfect cells according to the following table. The transfection is designed for 1 DNA amount combined with 4 different amounts of Lipofectamine® 2000. For each lipid reagent amount, the prepared mix is enough to have triplicates (96-well), duplicates (24-well), and single well (6-well) transfections, and account for pipetting variations.

Component	96-well	24-well	6-well
Adherent cells	1–4 x 10 <sup>4</sup>	0.5–2 x 10 <sup>5</sup>	0.25–1 x 10 <sup>6</sup>
Opti-MEM® Medium	25 µL x 4	50 µL x 4	150 µL x 4
Lipofectamine® 2000 Reagent	1, 1.5, 2, 2.5 µL	2, 3, 4, 5 µL	6, 9, 12, 15 µL
Opti-MEM® Medium	125 µL	250 µL	700 µL
DNA (0.5–5 µg/µL)	2.5 µg	5 µg	14 µg
Diluted DNA	25 µL	50 µL	150 µL
Diluted Lipofectamine® 2000 Reagent	25 µL	50 µL	150 µL
Incubate for 5 minutes at room temperature			
DNA-reagent complex/well	10 µL	50 µL	250 µL
Incubate cells for 1–3 days at 37°C			

The following table shows the amounts of DNA and Lipofectamine® 2000 Reagent per well used in each transfection reaction. For additional information on scaling your transfection reaction, see page 4.

Amount	96-well	24-well	6-well
DNA/well	100 ng	500 ng	2500 ng
Lipofectamine® 2000 Reagent/well	0.2–0.5 µL	1–2.5 µL	5–12.5 µL

## Scaling Up or Down Transfections

Use the following table to scale the volumes for your transfection experiment.

Culture Vessel	Multi- plication factor <sup>1</sup>	Shared reagents		DNA transfection		RNAi transfection	
		Vol. growth medium	Opti-MEM/ medium vol. for complex	DNA ( $\mu$ g)	Lipid reagent <sup>2</sup> ( $\mu$ L)	RNA (pmol)	Lipid reagent <sup>2</sup> ( $\mu$ L)
96-well	0.17	100 $\mu$ L	2 $\times$ 5 $\mu$ L	0.1	0.2–0.5	3	0.3
48-well	0.50	250 $\mu$ L	2 $\times$ 12.5 $\mu$ L	0.25	0.5–1.3	7.5	0.75
<b>24-well</b>	<b>1.00</b>	<b>500 <math>\mu</math>L</b>	<b>2 <math>\times</math> 25 <math>\mu</math>L</b>	<b>0.5</b>	<b>1–2.5</b>	<b>15</b>	<b>1.5</b>
12-well	2.00	1 mL	2 $\times$ 50 $\mu$ L	1	2–5	30	3
6-well	5.00	2 mL	2 $\times$ 100 $\mu$ L	2.5	5–12.5	75	7.5
60-mm	11.05	5 mL	2 $\times$ 250 $\mu$ L	5.5–11	11–28	166	17
10-cm	28.95	10 mL	2 $\times$ 500 $\mu$ L	14–28	29–73	434	43
T75	39.47	15 mL	2 $\times$ 750 $\mu$ L	20–40	39–100	592	59
T175	92.11	35 mL	2 $\times$ 1.75 mL	46–90	92–230	1382	138

<sup>1</sup>After determining the optimum reagent amount, use the multiplication factor to determine the reagent amount needed for your new plate format.

<sup>2</sup>Optimum amount needed is determined from the protocol (see pages 2–3).

## Co-Transfection of Plasmid DNA and siRNA

Transfect plasmid DNA and siRNA at the same time using Lipofectamine<sup>®</sup> 2000 Reagent by adding 30 pmol (~0.6  $\mu$ g) of siRNA per 1  $\mu$ g of DNA.

## mRNA Transfection

mRNA can be transfected in a 24-well plate by using Lipofectamine<sup>®</sup> 2000 Reagent by adding 0.5–1  $\mu$ g of mRNA per well.

### Limited Product Warranty

Life Technologies Corporation and/or its affiliate[s] warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at [www.lifetechnologies.com/terms](http://www.lifetechnologies.com/terms) and conditions. If you have any questions, please contact Life Technologies at [www.lifetechnologies.com/support](http://www.lifetechnologies.com/support).

**Disclaimer:** LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF.

©2012 Life Technologies Corporation. All rights reserved. The trademarks mentioned herein are the property of Life Technologies Corporation or their respective owners.

For support visit [www.lifetechnologies.com/support](http://www.lifetechnologies.com/support)

[www.lifetechnologies.com](http://www.lifetechnologies.com)

